

Serial number.....

Production year



SPÓŁDZIELNIA PRACY
46 Boremlowska Street
04-347 Warsaw/Poland
Phone (+48 22) 610 81 07 Service
Fax (+48 22) 610 55 36

10-04-2006

USER MANUAL

Catalogue number 20223es/ENG

LABORATORY CENTRIFUGE

MPW – 223 e
(10.223es – 5000 rpm)



Read it before use!

CE IVD

Contents

- 1. Application**
- 2. Technical data**
 - 2.1. Accessories
 - 2.1.1. Basic accessories
 - 2.1.2. Optional accessories
 - 2.2. Exploitation materials
- 3. Installation**
 - 3.1. Unpacking of the centrifuge
 - 3.2. Location
 - 3.3. Connection to mains
 - 3.4. Fuses
- 4. Description of the centrifuge**
 - 4.1. General description
- 5. Safe working conditions**
 - 5.1. Operating personnel
 - 5.2. Guarantee period and operation life
 - 5.3. Safekeeping period
 - 5.4. Hints on centrifuging
 - 5.5. Hazards and precautions
- 6. Operation of the centrifuge**
 - 6.1. Mounting of the rotor and accessories
 - 6.2. Construction and safety measures
 - 6.3. Drive
 - 6.4. Data setting and read-out
 - 6.5. Controls
 - 6.6. Safety devices
 - 6.6.1. Cover lock
 - 6.6.2. Rest state check
- 7. Description of the centrifuge operating elements**
 - 7.1. Control panel - Drawing No. 2
 - 7.2. Switching the centrifuge on
 - 7.2.1. Selection of the program
 - 7.2.2. Start of the program
 - 7.2.3. Emergency stop
 - 7.2.4. End of the centrifuging
 - 7.2.5. Programming
 - 7.3. Mathematical relations

- 7.3.1. RCF – relative centripetal force
- 7.3.2. Nomogram of relationship – rotational speed/centrifuging radius/RCF – Drawing No. 3
- 7.3.3. Maximum load

8. Cleaning, disinfection, maintenance

- 8.1. Cleaning of the centrifuge
- 8.2. Cleaning of the accessories
- 8.3. Lubrication
- 8.4. Glass cracking
- 8.5. Sterilization and disinfection of the rotating chamber and accessories

9. Emergency conditions – service

- 9.1. Troubleshooting

10. Safety work

- 10.1. Safety work inspection procedures
- 10.2. Inspection procedures carried out by the operator

11. Conditions of repairs

12. Manufacturer's data

13. Information about Distributor

1. Application.

The MPW-223e (the MPW-223es option) centrifuge is a table top laboratory centrifuge intended for *in vitro* diagnostic (IVD). Its construction ensures easy operation, safe work and wide range of applications in laboratories engaged in routine medical analyses, biochemical research works etc. It is intended for separation of mixtures, suspensions and systemic fluids into constituents of different densities under influence of the centrifugal force. This centrifuge is not biotight and therefore during centrifugation of preparations that require biotightness one has to use closed and sealed buckets and rotors. It is prohibited to centrifuge caustic, inflammable and explosive preparations in the centrifuge.

2. Technical data.

Manufacturer:

**“MPW MED, INSTRUMENTS”
SPÓŁDZIELNIA PRACY
46 Boremlowska Street, Warsaw/Poland**

Type:

MPW-223e Catalog No. 10223es

Mains: L1+N+PE V/Hz $\pm 10\%$	230 V 50/60 Hz, optionally 115 V 50/60 Hz
Power consumption	135 W
Class of protection against electric shock	basic
Interference level	PN-EN-55011
Noise level	58 dB
Rotational speed range	300÷5000 rpm
Maximum capacity	180 ml
Maximum acceleration RCF	3689 x g
Maximum kinetic energy	1536 Nm
Time range	1 ÷ 99 min
SHORT – short duration operation	

Physical data:

Depth	435 mm
Width	355 mm
Height	270 mm
Weight	13 kg

Centrifuge operation conditions:

PN-EN 61010-1 p. 1.4.1.

Environmental temperature	+2° ÷ +40° C
Relative humidity at ambient temperature	< 80%
Installation category	II PN-EN 61010-1
Degree of pollution	2 PN-EN 61010-1
Protection zone	300 mm

Statement of Conformity:

The following machine is in accordance with the regulations of the EU Directive 98/79/EC and with the harmonized standards PN-EN 61010-1 and PN-EN 61010-2-020.

2.1. Accessories.

2.1.1. Basic accessories (enclosed with every centrifuge).

- 17099T	spanner for the rotor	pcs. 1
- 17162	key for the rotor	“ 1
- 17142	complete clamp	“ 1
- 17861	fuses WTA-T 4 A 250 V	“ 2
- 17866	power cord 230 V	“ 1
- 17867	power cord 115 V (optionally)	“ 1
- 20223es /ENG	Operating Instruction	“ 1

2.1.2. Optional accessories.

Depending on costumer's need the MPW-223e centrifuge can be provided with accessories specified below:

Indeks Cat. No	Nazwa Name MPW-223es	Probówka MPW MPW Tube
11237C/A	Wirnik kątowy 12 x 15/10ml, komplet z pojemnikami 13080 (O 17x100/120mm) (kąąt 30°) Angle rotor 12 x 15ml, complete with buckets 13080 (O 17x100/120mm) (angle 30°) (max RPM: 5000 max RCF: 3326 x g R max: 11,9 cm)	15048,15050, 15053,15118
14082	Wkładka redukcyjna (O 13,3mm) Round carrier (O 13,3mm)	15119
14083	Wkładka redukcyjna (O 13,3mm) **** DO WYCZERPANIA ZAPASÓW **** Round carrier (O 13,3mm) **** SELL OUT ****	15119
11237C/B	Wirnik kątowy 12 x 10ml, komplet z pojemnikami 13081 (O 17x70/85mm) (kąąt 30°) Angle rotor 12 x 10ml, complete with buckets 13081 (O 17x70/85mm) (angle 30°) (max RPM: 5000 max RCF: 3326 x g R max: 11,9 cm)	15053
14082	Wkładka redukcyjna (O 13,3mm) Round carrier (O 13,3mm)	15054,15120, 15419
14083	Wkładka redukcyjna (O 13,3mm) **** DO WYCZERPANIA ZAPASÓW **** Round carrier (O 13,3mm) **** SELL OUT ****	15054,15120, 15419
11257C	Wirnik kątowy 4 x 30/25ml, komplet z pojemnikami 13329 (kąąt 45°) Angle rotor 4 x 30/25ml, complete with buckets 13329 (angle 45°) (max RPM: 5000 max RCF: 3326 x g R max: 11,9 cm)	15055, 15056, 15424
14255	Wkładka redukcyjna na probówkę 7ml (O 13x100mm) Round carrier for 7ml tube (O 13/100mm)	15054, 15119
14256	Wkładka redukcyjna na probówkę 15/10ml (O 17x120mm) Round carrier for 15/10ml tube (O 17/120mm)	15046, 15048, 15053, 15118
11487C/A	Wirnik kątowy 8 x 15/10ml, komplet z pojemnikami 13080 (O 17x100/120mm) (kąąt 45°) Angle rotor 8 x 15/10ml, complete with buckets 13080 (O 17x100/120mm) (angle 45°) (max RPM: 5000 max RCF: 3214 x g R max: 11,5 cm)	15048,15050, 15053,15118
14082	Wkładka redukcyjna (O 13,3mm) Round carrier (O 13,3mm)	15119
14083	Wkładka redukcyjna (O 13,3mm) **** DO WYCZERPANIA ZAPASÓW **** Round carrier (O 13,3mm) **** SELL OUT ****	15119
11487C/B	Wirnik kątowy 8 x 10ml, komplet z pojemnikami 13081 (O 17x70/85mm) Angle rotor 8 x 10ml, complete with buckets 13081 (O 17x70/85mm) (angle 45°) (max RPM: 5000 max RCF: 3214 x g R max: 11,5 cm)	15053
14082	Wkładka redukcyjna (O 13,3mm) Round carrier (O 13,3mm)	15054,15120, 15419

14083	Wkładka redukcyjna (O 13,3mm) **** DO WYCZERPANIA ZAPASÓW **** Round carrier (O 13,3mm) **** SELL OUT ****	15054,15120, 15419
11720C	Wirnik kątowy 4 x 60ml na probówki Orthokine®vet, komplet z pojemnikami 13721 Angle rotor 4 x 60ml for Orthokine®vet tubes, complete with buckets 13721 (max RPM: 5000 max RCF: 3633 x g R max: 13 cm)	
11725C	Wirnik kątowy 2 x 50ml + 4 x 20ml, komplet z 2 pojemnikami 13726 na strzykawki BD PERFUSION 50ml SYRINGE® oraz z 4 pojemnikami 13727 na strzykawki KD-JECT III 20ml® (kąt 45°) Angle rotor 2 x 50ml + 4 x 20ml, complete with two 13726 buckets for BD PERFUSION 50ml SYRINGE® and with four 13727 buckets for KD-JECT III 20ml® syringe (angle 45°) (max RPM: 5000 max RCF: 3689 x g R max: 13,2 cm)	

Indeks Article symbol	Nazwa Name Probówki Test tubes
15048	Probówka z polipropylenu 15ml Nalgene® (O 16x113mm) Polypropylene tube 15ml Nalgene® (O 16x113mm)
15050	Probówka z polipropylenu 15ml z dnem stożkowym (O 17/21x120mm) Polypropylene tube 15ml with conical bottom
15053	Probówka z polipropylenu 10ml z pokrywką (O 16x100mm) Polypropylene tube 10ml with cap (O 16x100mm)
15054	Probówka z polipropylenu 6ml z pokrywką (O 11,7/13,5x95mm) Polypropylene tube 6ml with cap (O 11,7/13,5x95mm)
15055	Probówka z polipropylenu 30ml z pokrywką (O 24,9x103mm) Polypropylene tube 30ml with cap (O 24,9x103mm)
15056	Probówka z poliwęglanu 30ml Nalgene® z pokrywką (O 25,5x94mm) Polycarbonate tube 30ml Nalgene® with cap (O 25,5x94mm)
15118	Probówka szklana 10ml (O 16x100mm) Glass tube 10ml (O 16x100mm)
15119	Probówka szklana 7ml (O 12x100mm) Glass tube 7ml (O 12x100mm)
15120	Probówka szklana 5ml (O 12x75mm) Glass tube 5ml (O 12x75mm)
15419	Probówka z polipropylenu 5ml z korkiem (O 12x85mm) Polypropylene tube 5ml (O 12x85mm) with cap
15424	Probówka z polipropylenu 30ml z pokrywką (O 25,5x94mm) Polypropylene tube 30ml with cap (O 25,5x94mm)

2.2. Exploitation materials.

For centrifuge operation one should use only original company brand buckets comprised in the specification of accessories as well as test-tubes for centrifuges of proper diameter, length and strength. Use of test-tubes of other makes shall be agreed upon with manufacturer of the centrifuge. For cleaning and disinfecting one should to use agents generally used in the health service, such as e.g. Aerodesina-2000, Lysoformin 3000, Malseptol, Malsept SF, Sanepidex, Cutasept F.

3. Installation.

3.1. Unpacking of the centrifuge.

Open the package. Take out the cardboard box containing the accessories. Take out the centrifuge from the package. Keep the package and packing materials at hand for possible transport at a later date.

3.2. Location.

Almost all energy being supplied to the centrifuge is transformed into heat and then emitted to the environment. This is the reason why proper ventilation is essential. Ventilation ducts situated in the centrifuge have to be fully efficient. Moreover the centrifuge shall not be located near the radiators and shall not be subjected to direct sunlight. The table for the centrifuge shall be stable and shall have flat levelled table top. Safety zone has to be established around the centrifuge with the minimum radius of 30 cm. Under normal operating conditions ambient temperature shall not drop below 15° C or raise above 35° C. In the case of changing location from a cold to warm one there will occur condensation of water inside the centrifuge. It is important then that sufficient time shall be allowed for drying the centrifuge prior to restarting of the centrifuge (minimum 4 hours).

3.3. Connection to mains.

Supply voltage given on the rating plate has to be consistent with local supply voltage.

MPW Med. instruments laboratory centrifuges are I safety class devices and are provided with the three-core cable of 2.5÷3.2 m length with the plug resistant to dynamic loadings. Mains socket shall be provided with the safety pin. Neutral grounding of mains socket safety pin has to be verified by authorized services. This verification has to be carried out each time when mains socket is being replaced. It is recommended to install emergency cut-out that shall be installed far from the centrifuge, near the exit door from the room or outside the room. Supply voltage 230 V 50/60 Hz, optionally 115 V 50/60 Hz.

3.4. Fuses.

The centrifuge has standard protection with the WTA-T 4A 250 V fuse situated in the plug-in socket and master switch unit at back wall of the centrifuge.

4. Description of the centrifuge.

4.1. General description.

New generation of MPW Med. instruments laboratory centrifuges is provided with modern microprocessor control systems, very durable and quiet asynchronous brushless motors and accessories consistent with modern requirements.

5. Safe working conditions.

5.1. Operating personnel.

MPW-223e Laboratory Centrifuge can be operated by laboratory personnel after getting acquainted with the Instruction Manual.

Instruction Manual shall be always near the centrifuge.

Instruction Manual must be constantly at hand!!

5.2. Guarantee period and operation life.

Guarantee period for MPW-223e centrifuge amounts to minimum 24 months.

Principles are specified in guarantee certificate. The service life of the centrifuge specified by the manufacturer amounts to 10 years.

After termination of guarantee period it is necessary to carry out annual technical inspections of the centrifuge made by authorized service of manufacturer.

The manufacturer reserves the right to make modifications at produced goods.

5.3. Safekeeping period.

Maximum period of storage of not used centrifuge amounts to 1 year. After this period one should ask authorized service to carry out an inspection of the device.

5.4. Hints on centrifuging.

1. Set the centrifuge in horizontal position on rigid base.
2. Ensure safe location.
3. Ensure free space around the centrifuge, within at least 30 cm.
4. Ensure sufficient ventilation.
5. Fix firmly the rotor on the motor shaft.

6. **Avoid unbalance.**
7. **Load opposite buckets with the same accessories.**
8. **Centrifugation of the test tubes of different dimensions.**
 In principle it is possible to centrifuge test tubes of different dimensions, however it is absolutely necessary in such cases that opposite round carriers have to be the same.
 The test tubes shall be not only inserted symmetrically but round carriers and their hangers shall be equally loaded. It is not allowed to operate centrifuge with asymmetric loads applied to rotors and buckets.
9. **Load all places in rotors.**
10. **Fill test tubes outside the centrifuge.**
11. **Glass tubes shall be test tubes intended for centrifuges of proper strength enabling centrifugation with acceleration up to 5000 x g.**
12. **Fill in the test tubes with the medium of the same weight, in order to protect the centrifuge against unbalance.**
13. **Lubricate the rotor journal pins.**
14. **Use only accessories kept in good condition.**
15. **Protect equipment against corrosion using accurate preventive maintenance.**
16. **Infectious materials could be processed in closed buckets only.**
17. **It is prohibited to centrifuge explosive and inflammable materials.**
18. **It is prohibited to centrifuge substances prone to reacting in result of supplying high energy during centrifugation.**

5.5. Hazards and precautions.

1. **Prior to starting the trial of switching the centrifuge on, one shall read exactly all sections of this instruction in order to ensure smooth run of operation, avoiding damages of this device or its accessories.**
2. **Centrifuge can be operated by laboratory personnel getting acquainted with the Instruction Manual.**
3. **Centrifuge must not be transported with the rotor mounted on the motor shaft.**
4. **One must use original rotors, test-tubes and spare parts only.**
5. **In case of faulty operation of the centrifuge one shall ask for assistance of service of MPW Med. instruments Company or its accredited representatives.**
6. **It is prohibited to switch the centrifuge on if it is not installed properly or rotor is not fitted correctly.**
7. **The centrifuge must not be operated in places where explosion hazard appears as it is not of explosion-proof make.**
8. **It is prohibited to centrifugate materials which could generate inflammable or explosive mixtures when in contact with air.**
9. **It is prohibited to subject to centrifugation toxic or infectious materials without taking proper safety measures (work in properly adapted rooms, personal safety equipment). Proper disinfection procedures have to be carried out when dangerous substances contaminated the centrifuge or its accessories.**
10. **One must not open the cover manually – in emergency procedure, when rotor is still turning.**
11. **One must not exceed limit load set by the manufacturer.**
 Rotors are intended for fluids of average homogeneous density equal to 1.2 g/cm³ or smaller when centrifugation is carried out at maximum speed. When fluids of higher density shall be used, then it is necessary to limit the speed (see point 7.3.3 “Maximum load”).
12. **One must not to use the rotors, buckets and round carriers with symptoms of corrosion or other mechanical defects.**
13. **One must not subject to centrifugation substances of high corrosion aggressiveness, which could cause material impairment and lower mechanical properties of rotors, buckets and round carriers.**
14. **One must not use rotors and accessories not admitted by the manufacturer.**
 It is permitted however to use commercial glass and plastic test tubes intended by manufacturer of these test tubes for centrifugation in laboratory centrifuges. It is distinct warning against using

accessories not specified in the Instruction Manual. Cracking of test tubes not intended for centrifugation can result in dangerous unbalance.

15. One must not carry out centrifugation with the rotors with taken off or not tight driven caps.
16. One must not lift or shift the centrifuge during operation and rest on it.
17. One must not stay in the safety zone within 30 cm distance around the centrifuge neither leave any things, e.g. glass vessels, within this zone.
18. It is prohibited to put any things on the centrifuge.

6. Operation of the centrifuge.

6.1. Mounting of the rotor and accessories.

1. Connect the centrifuge to mains (master switch at back wall of the centrifuge).
2. Open the cover of the centrifuge pushing the pushbutton COVER. Prior to putting the rotor in one has to check if rotating chamber is free of impurities, e.g. such as dust, glass splinters, residues of fluids that must be taken away.
3. One shall release with special spanner clamp on the motor shaft and fit the rotor on the motor shaft driving it home on the cone.
4. Screw-in the bolt for fixing the rotor (clockwise) and screw it tightly home with the supplied spanner.
5. In the case of rotors designed with the cover they must not be used without it. Rotor covers must be closed exactly. Rotor covers ensure smaller drags of the rotors, proper setting of the test-tubes and airtight sealing.
6. One should use only buckets intended for selected types of the rotor - see p. 2.1. "Accessories".
7. Fill test tubes outside the centrifuge.
8. Put on or screw the caps on vessels and rotors (when they have such).
9. In the case of centrifuging in angle rotor, test tubes (buckets) have to be filled properly in order to avoid overflows.
10. **CAUTION:** *Centrifuge will tolerate small weight differences occuring during loading of rotors. However it is recommended to equalize vessels loads as much as possible in order to ensure minimal vibrations during operation.*
11. Maximum unbalance amounts up to 15 g.
12. For replacement of the rotor one shall release clamping by several turns of the bolt and then using both hands grab the rotor at opposite sides taking it away from drive shaft by pulling it up.

6.2. Construction and safety measures.

The centrifuge has rigid self-supporting structure. Housing is made of plastic while front is made of steel sheet. Cover is fixed on steel axles of hinges and from the front is locked with electromagnetic lock blocking possible opening during centrifugation. Bowl forming the rotation chamber is made of acid resistant steel sheet.

6.3. Drive.

Drive constitutes brushless induction motor of low noise level, free of carbon brushes. This solution eliminated the danger of contamination the preparations with carbon dust.

6.4. Data input and output.

Data setting and read-out system forms hermetically closed keyboard with distinctly accessible operation points. Easily readable display signalling individual performed operations facilitates to operator programming of condition of the centrifuge.

Operation of the centrifuge is simple and self-evident.

6.5. Controls.

Microprocessor control system being used in the centrifuge ensures the following possibilities of setting of parameters of operation:

- stepwise rotational speed selection within the range from 300 up to 5000 rpm every 100 rpm with calculating and RCF readout,
- setting of the centrifugation time within the range from 1 up to 99 minutes.

6.6. Safety devices.

Apart from the above described passive devices and safety measures there exist as well active devices and elements as follows:

6.6.1. Cover lock.

The centrifuge can be started only with properly closed cover. The cover can be opened only after stopping the rotor. In the case of emergency opening of the cover during operation the centrifuge will be immediately switched-off and the rotor will be braked until stopping completely. With opened cover, the drive is completely disconnected from power which makes it impossible to start the centrifuge.

6.6.2. Rest state check.

Opening of the centrifuge cover is possible only with the rotor in the state of rest. This state is being checked by the microprocessor which recognizes and signals the rest state prior to opening the cover with letter **S** (Stop).




7. Description of the centrifuge operating elements.

Power switching ON/OFF is carried out with master switch situated on back of the centrifuge. All settings on the centrifuge are done by means of the control panel. Panel comprises control pushbuttons, displays and signalling LEDS.


7.1. Control panel.


Control panel placed on front casing wall serves for controlling centrifuge operation. Control panel comprises the following elements:




- | | |
|----------------------------|--|
| 1. Upper display field | S: 4 digits (rpm) RCF 4 digits (x g) |
| 2. Lower display field | T: 4 digits (m/s) S (STOP), O (opened cover), U (unbalance) |
| 3. Error signalling | |
| 4. Rotor status signalling | STATUS COVER blinking - rotor rotates,
not illuminated - rotor does not rotate |


- | | |
|-----------------|---|
| 5. Function key |  |
| 6. Function key |  |
| 7. Function key |  |

- ▶ Buzzer serves for signalling function recording and determination of the centrifuge status.

- ▶  key serves for starting centrifugation program with parameters presented on display,

- ▶  key serves for:
 - interrupting centrifugation program in any program phase and braking the rotor,
 - saving of preset SPEED and TIME centrifugation parameters,

- ▶  key serves for opening of the cover,
- ▶  serve for speed programming,
- ▶  serve for time programming,

- ▶  key serves for short duration operation.

7.2. Switching the centrifuge on.

After switching power ON the control system calls recently implemented program and displays in relevant fields rotational speed, duration of centrifugation and cover opening status. Provided that rotor in the centrifuge is stopped, it is possible to open the cover by means of COVER key.

Stopped rotor status is displayed as a **S** letter symbol in the display field. When this symbol is not already displayed, then one must wait till this rotor stops and the above mentioned symbol appears.

7.2.1. Selection of the program.

Control panel can save 1 program preset by the user. Program acceptance consists in pushing STOP key.

7.2.2. Start of the program.

After acceptance of the program and checking if rotor was mounted, centrifuge can be started with pushing START key, provided that cover is closed.

7.2.3. Emergency stop.

At any time during centrifuging it is possible to interrupt the process and stop the rotor quickly with single pushing of STOP key.

7.2.4. End of the centrifuging.

After termination of time of centrifugation preset in the program, braking follows. At the end of deceleration the rotational speed drops at slower rate in order to ensure soft settling of rotor carriers. Stopping is followed by buzzer signal and is displayed by **S** letter symbol. After pushing COVER pushbutton the cover opens and **O** symbol is displayed.

7.2.5. Programming.

Programming mode is activated with pushing SPEED and TIME (+) (-) keys after selection of parameters of the program which one would like to save or change. Acceptance of preset parameters is done by pushing STOP key. One can save one program only.

7.3. Mathematical relations.

7.3.1. RCF - relative centripetal force.

RCF acceleration is the acceleration generated by the rotor rotary motion acting upon tested product and it can be calculated according to the formula:

$$RCF = 11.18 * r * (n/1000)^2$$

$$\begin{array}{lll} \text{where} & RCF & [x \text{ g}] \\ r & [cm] \\ n & [rpm] \end{array}$$

Depending on the distance of particles of the tested product from the axis of rotation one can find from above formula minimum RCF, average RCF or maximum RCF. On the basis of preset RCF value and given radius of the bottom in the bucket one can calculate from the formula the rotational speed to be set in the program of centrifuging. Selection of the time of sedimentation and the RCF value shall be carried out experimentally for a given product.

Once every 100 rpm electronic circuit automatically calculates and displays averaged RCF value.

7.3.2. Nomogram of relationship – rotational speed/centrifuging radius/RCF – Drawing No. 3

7.3.3. Maximum load.

In order to avoid overloading of the rotor one shall observe maximum load which is recorded on every rotor. Maximum permissible load is reached when all test-tubes are filled with the fluid with 1.2 g/cm³ density.

If density of the centrifugated liquid is higher than 1.2 g/cm³, then test-tubes could be filled only partially or one shall limit operation speed of the centrifuge that is being calculated from the formula:

$$n_{perm} = n_{max} * \sqrt{\frac{1,2}{\gamma}}$$

$$\gamma = \text{specific gravity} \left[\frac{G}{cm^3} \right]$$

n_{max} [maximum rotational speed – rpm]

8. Cleaning, disinfection, maintenance.

CAUTION!!! Use safety gloves for operations specified below.

8.1. Cleaning of the centrifuge.

Prior to start of cleaning and disinfection of the centrifuge one shall put the safety gloves on. For cleaning shall be used water with soap or other water soluble mild detergent. One should avoid corrosion inducing substances and aggressive substances. It is prohibited to use alkaline solutions, inflammable solvents or agents containing abrasive particles.

8.2. Cleaning of the accessories.

In order to ensure safety operation one shall in regular way carry out periodical maintenance of the accessories. Manufactured rotors, buckets and round carriers have to withstand high stresses originated from centrifugal force. Chemical reactions as well as corrosion (combination of variable pressure and chemical reactions) can cause corrosion or destruction of metals. Hard to observe surface cracks increase gradually and weaken material without visible symptoms. In the case of observation of surface damage, crevice or other change, as well the corrosion, given part (rotor, bucket, etc.) shall be immediately replaced. In order to prevent corrosion one has to clean regularly the rotor together with the fastening bolt, buckets and round carriers. Cleaning of the accessories shall be carried out outside of the centrifuge once every week or still better after each use. Then those parts shall be dried using soft fabric or in the chamber drier at ca. 50° C. Especially prone to the corrosion are parts made of aluminium. For cleaning them one should use very neutral agent of pH value within 6÷8 range. It is forbidden to use alkaline agents of pH above 8. In this way substantially is increased useful service life and diminished susceptibility to corrosion. Accurate maintenance also increases service life and protects against premature rotor failures. Corrosion and damages resulting from insufficient maintenance could not be object of claims lodged against the manufacturer.

8.3. Lubrication.

The rotor pins shall be always lubricated with acid-free vaseline. In this way is ensured uniform deflection of the buckets and quiet centrifuge operation.

8.4. Glass cracking.

In the case of glass cracking one shall wear safety gloves and all debris shall be accurately removed. Rubber inserts shall be exactly cleaned or possibly replaced. Otherwise one has to take into account the following possibilities:

- Glass particles left in the rubber cushion (pad) will cause once more glass cracking.
- Glass particles left in containers make impossible uniform deflecting of the buckets and round carriers resulting in unbalance.
- Glass particles left in the rotor chamber cause metal abrasion because of strong air circulation. This dust will not only contaminate the centrifuge chamber, rotor, buckets, carriers and centrifuged material but will cause as well damages of surfaces of the accessories, rotors and the centrifuge chamber. For complete removal of glass particles and metal dust from the rotor chamber it is recommended to place on the bowl strip of vaseline (from the top down to bottom). Then rotor shall operate for several minutes at moderate speed. Glass and metal particles will collect on lubricated area and could be easily removed with the piece of cloth together with the grease. This operation can be repeated in case of need.

8.5. Sterilization and disinfection of the rotating chamber and accessories.

One can use all standard disinfectants. The centrifuges and accessories are constructed from various materials and one should take into account possible variety of materials. During sterilization by means of steam one should consider temperature resistance of individual materials. We would like to add that for centrifuging for instance infectious materials it is necessary to use hermetically closed buckets in order to protect their migration into the centrifuge. Buckets and rotors can be sterilized in autoclave with temperature 121°–124° C and pressure 215 kPa during 15 min. Disinfection is carried out with disinfectants generally used in the Health Service (see paragraph 2.2.).

User is responsible for proper disinfection of the centrifuge, if some dangerous material was spilled inside or outside the centrifuge. During above mentioned works one must wear safety gloves.

9. Emergency conditions – service.

9.1. Troubleshooting.

Majority of faults could be cancelled by switching the centrifuge OFF and then ON. After switching the centrifuge ON shall be displayed parameters of the recently implemented program and buzzer signals consisting of four successive tones. In the case of short-duration power failure the centrifuge terminates cycle.

Please find below the most frequent faults and their repair methods.

1. Lack of display and check buzzer:	Remedies:
<i>Is mains socket live?</i>	<i>Check mains socket fuse.</i>
<i>Is supply cable plugged into mains?</i>	<i>Plug correctly supply cable.</i>
<i>Is input fuse good?</i>	<i>Replace input fuse (rated data on rating plate).</i>
<i>Is master switch switched ON?</i>	<i>Switch ON power supply.</i>
<i>The above was checked and still there is no display active and no check buzzer sound.</i>	<i>Call service.</i>
2. Centrifuge does not start	Remedies
<i>START key pushing does not generate reaction or single tone only</i>	
<i>Rotor stopping symbol S is not displayed yet</i>	<i>Wait till rotor stops and the rotor stopping symbol is displayed</i>
<i>Cover opening symbol O is displayed</i>	<i>Close cover. S symbol that means stop should be displayed.</i>
<i>LED STATUS diode is blinking:</i>	<i>Centrifugation cycle in progress, push STOP key or wait till cycle ends.</i>
<i>Indications show a cycle in progress but the motor does not start</i>	<i>Switch power supply OFF/ON. If fault still persists then call service.</i>
3. Programming function not active	Remedies
<i>It is impossible to record parameter values to memory, last recorded program can not be recalled. Disturbances on displays possible too.</i>	<i>Call service.</i>
4. Centrifuge starts but does not accelerate	Remedies
<i>E symbol displayed after stopping. Drive overload</i>	<i>Wait for 15 minutes and switch again after opening and closing the cover.</i>
5. One can not open the cover	Remedies
<i>Rotor stopping S symbol not displayed yet, after pushing cover opening key single tone is audible</i>	<i>Rotor is still rotating. Wait for stopping of the rotor and displaying of the S symbol.</i>
<i>Nothing is displayed</i>	<i>Check the centrifuge power supply</i>
<i>Rotor stopping S symbol is displayed, but cover cannot be opened</i>	<i>Call service</i>

Emergency cover release

In the possible case of failure, such e.g. as power failure there exists possibility of manual opening of the cover. On the right side of the housing there is a small opening in which one shall put $\phi 2\text{mm}$ bar or key and push it, and the cover will be opened by itself.

CAUTION! Cover can be released and opened only when rotor is in the rest state.

10. Safety work.

10.1. Safety work inspection procedures.

From the point of view of operational safety the centrifuge has to be subjected to inspection carried out by authorized service engineer or especially trained experts at least once every year in the state of operational readiness. The reason for more frequent inspection could be for instance more frequent unbalance cases or corrosion inducing environment. Results of inspections, repairs and tests have to be recorded and kept on file. Operating Instruction shall be stored in the centrifuge use place.

10.2. Inspection procedures carried out by the operator.

Operator has to pay special attention to the fact that the centrifuge parts important because of safety reasons are not damaged.

This remark is specifically important for:

1. Motor suspension
2. Motor shaft concentricity
3. Centrifuge accessories and especially structural changes, corrosion, preliminary cracks, abrasion of metal parts
4. Screw joints
5. Inspection of the rotor assembly
6. Inspection of protective joints.

11. Conditions of repairs.

Manufacturer grants to the Buyer a guarantee on conditions specified in the Guarantee Certificate. Buyer forfeits the right to guarantee repair when using the device inconsistently with the Operating Instruction provisions, when damage resulted from the User's fault. Repairs should be carried out in authorized service workshops granted with the MPW Certificate. The centrifuge shall be sent to repair after decontaminating disinfection.

Information about authorized service workshops could be obtained from the Manufacturer, i.e. "MPW MED. INSTRUMENTS" in Warsaw, 46 Boremlowska Street.

Phone (+ 48 22) 610 81 07 or Exchange (+48 22) 610 50 14.

12. Manufacturer's data.

"MPW MED. INSTRUMENTS" 04-347 Warsaw, 46 Boremlowska Street

Phone: (+48 22) 610 56 67

Exchange: (+48 22) 610 50 14

Fax: (+48 22) 610 55 26


Internet: <http://www.mpw.pl>

E-mail: mpw@mpw.pl

13. Information about Distributor.

YOUR DISTRIBUTOR:

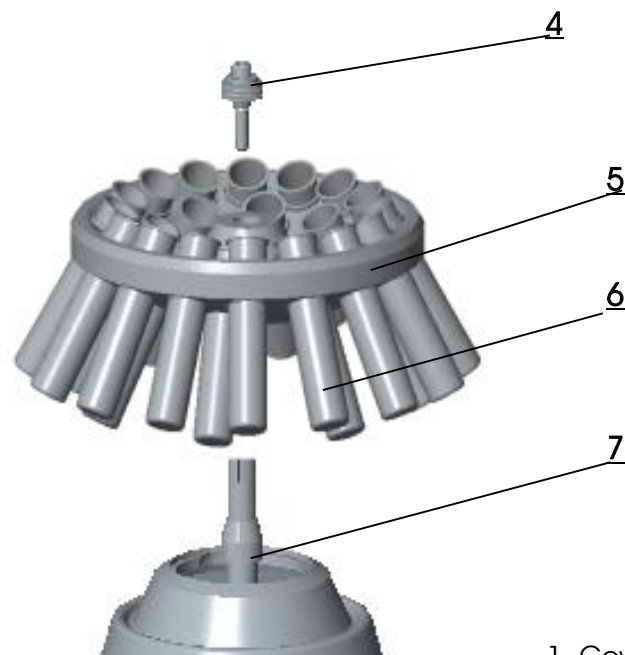


Drawing No. 1. General view.

1. Cover of the centrifuge
2. Emergency cover release
3. Control panel
4. Clamping bolt
5. Rotor
6. Bucket
7. Motor shaft

Drawing No. 2. Control panel.

Drawing No. 3. Nomogram.

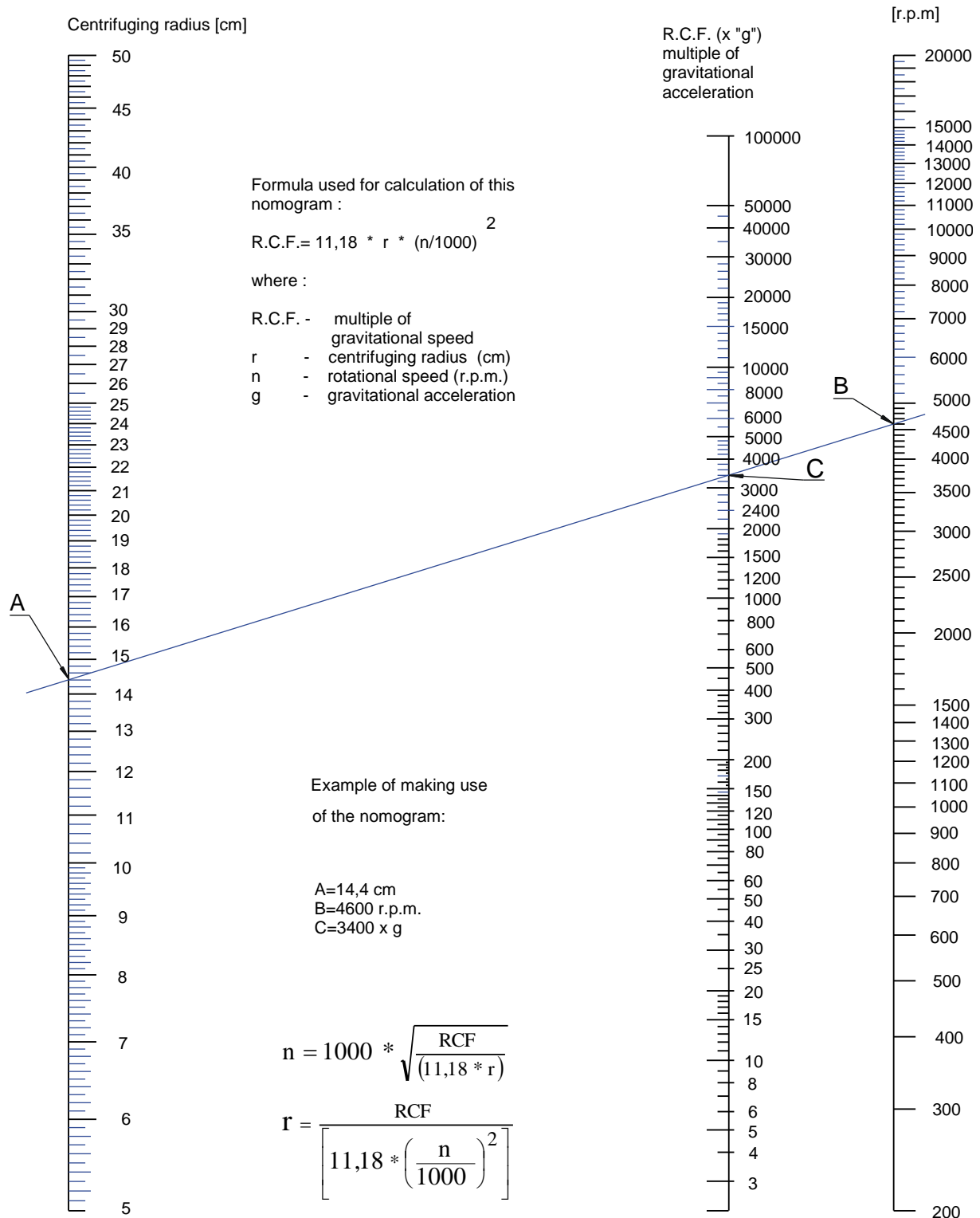

DECLARATION OF CONFORMITY

Product

Laboratory centrifuge

Model

MPW-223es

Product classification on the basis of
the Directive 98/79/EC

Non classified to list A or B and not for
self-testing

Product complies with the requirements:

• **Directive 98/79/EC (IVD), including the requirements of harmonised standards:**

PN-EN ISO 13485:2012

PN-EN ISO 18113-3:2011

PN-EN ISO 13485:2012/AC:2013-03

PN-EN 61010-2-101:2005

PN-EN 13612:2006

PN-EN 61326-2-6:2006

PN-EN ISO 14971:2012

PN-EN ISO 62366:2008

• **selected harmonized standards of Directive 2006/95/EC (LVD):**

PN-EN 61010-1:2011

PN-EN 61010-2-020:2008

• **Directive 2004/108/WE (EMC)**

• **standard PN-EN ISO 15223-1:2012**

CZŁONEK ZARZĄDU PREZES ZARZĄDU

Wojciech Nojszewski

mgr Hanna Matyńska

**„MPW MED. INSTRUMENTS”
SPÓŁDZIELNIA PRACY
w Warszawie**

**„MPW MED. INSTRUMENTS”
SPÓŁDZIELNIA PRACY**

Warsaw, 46 Boremlowska Street
Quality policy in line with ISO 9001:2008
Certifying authority



Warsaw, 21.05.2015

no 10.223es.03

DECLARATION OF DECONTAMINATION

In order to protect our employees please fill out the declaration of decontamination completely before sending centrifuge to the manufacturer (repair).

1. **Device**

– type:

– serial No.:

2. **Description of decontamination**

(see user manual)

.....

.....

.....

.....

3. **Decontamination carried out by:**

– name:

4. **Date and signature**

.....

DECLARATION OF DECONTAMINATION

In order to protect our employees please fill out the declaration of decontamination completely before sending centrifuge to the manufacturer (return).

5. **Device**

– type:

– serial No.:

6. **Description of decontamination**

(see user manual)

.....

.....

.....

.....

7. **Decontamination carried out by:**

– name:

8. **Date and signature**

.....